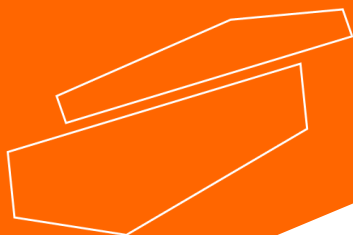
EVA180 Rotary evaporator

Please read the User Manual carefully before use, and follow all operating and safety instructions!

user manual

english / español / français / italiano

User Manual

LBX EVA180 Rotary evaporator

Preface

Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

Service

If help is needed, you can always contact your dealer or Labbox via www.labbox.com (declare an incidence)

Please, provide the customer service representative with the following information:

- Serial number (on the back side)
- Description of the problem
- Your contact information

Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

1. Safety instructions

	Connect the equipment to a grounded power source to ensure the safety of the instrument and the experiment; connect power as per the equipment requirements.
	The use of this equipment in flammable, explosive, toxic or highly corrosive experiments is prohibited.
	Place the equipment on a horizontal, flat and stable table, creating a free space of 30 cm per side, allowing you to work comfortably. Make sure the surface is adequate for the weight of the equipment.
	The equipment must be used by qualified personnel who are familiar with the equipment and its operation through the user manual.
	Do not place the equipment in areas close to heat sources.
	While the machine is in operation, do not touch the heating surface to avoid burns.
	During operation, hazardous material such as flammable liquids or pathological material must be away from this area.
	Read the instruction manual before using the equipment and keep it on a place on which it can be accessed by anyone.
	Use the appropriate safety elements, such as eye protection, gloves, according to the information provided by the reagent's manufacturer.
	Pay special attention to splashes and protect yourself adequately.

-During the use of the equipment, use personal protective measures to avoid risks of possible damage such as:

- Splash burns and evaporation of liquids
- Poisoning by the emission of toxic or combustible gases
- Cuts or physical damage due to the use of fragile material such as glass.

-Beware of the dangers caused by:

- Flammable materials or media with a low boiling temperature
- Overfilling of the container
- Unsafe containers

- Make sure that the glass assembly has no stress due to the risk of breaking from causes as:
 - Mechanical hazards from the outside
 - Excess temperature in the equipment itself
- The temperature must be always set, at least, 25°C below the ignition temperature of the materials used
- Check equipment and accessories before each use and do not use components in poor condition
- Safe operation of the equipment is only guaranteed with the accessories described in chapter "Packing list". The accessories must be firmly attached to the equipment and cannot come loose
- Always turn off the power before installing accessories
- Do not cover the ventilation outlets of the unit to ensure adequate cooling of the rotary evaporator
- Adapt the quantity and type of solvent to the size of the distillation equipment
 - Please know that high concentrations of gases, steam or particular matters can escape through the cooler. Avoid this risk by using, for example, cold water traps, gas wash bottles or an effective extraction system, such as showcases
- The glassware is designed for operations with maximum vacuum of 10 mbar. The glass must be cooled down before its removal and reuse
 - Avoid formation of peroxide. Organic peroxides can accumulate in the distillation and generate exhaust residues that explode easily when decomposed
 - Never heat the evaporation flask in the bath without first actuating the rotary motion. In this way, you will avoid possible losses of your sample
 - Pay attention to the temperature setting; never leave the equipment unattended during operation and with the heating function on
 - Do not handle the rotary evaporator when the rotating piston is working. It is recommended to gradually increase and/or decrease the rotation speed if situations such as an irregular operation occur
 - The equipment can only be disconnected from the power outlet by pulling the plug base, not the cord
 - The voltage indicated on the equipment must correspond to the electrical network used
 - Make sure that the main power cord is not in contact with the heating surface. Do not cover the equipment
 - Keep the equipment away from high magnetic fields

2. Rules

This equipment is specially designed for distillations for educational, laboratory or industrial purposes. This apparatus is not suitable for use at home or in environments that could be hazardous to the user or the equipment.

With the accessories recommended by the manufacturer, this equipment is capable of the following actions:

- Rapid distillation of liquids
- Evaporation of solutions and suspensions
- Crystallization, synthesis or cleaning of chemicals
- Recycle solvents

3. Inspection

3.1 Reception

Carefully unpack the instrument and check that the equipment and/or accessories have arrived undamaged. If necessary, contact the supplier who supplied the equipment for technical assistance.

Note:

- If the equipment is damaged, it must not be connected to an electrical source.
- Ensure that the device is unlocked before turning it on.

3.2 Packing list

The package includes the following items:

Item	Quantity
Main Unit	1
Heating bath	1
Condenser	1
Distillation flask (29/32)	1
Collector flask (35/20)	1
Support rod	1
Power cable	1
User manual	1

Table 1

4. Control elements

4.1 Parts and controls

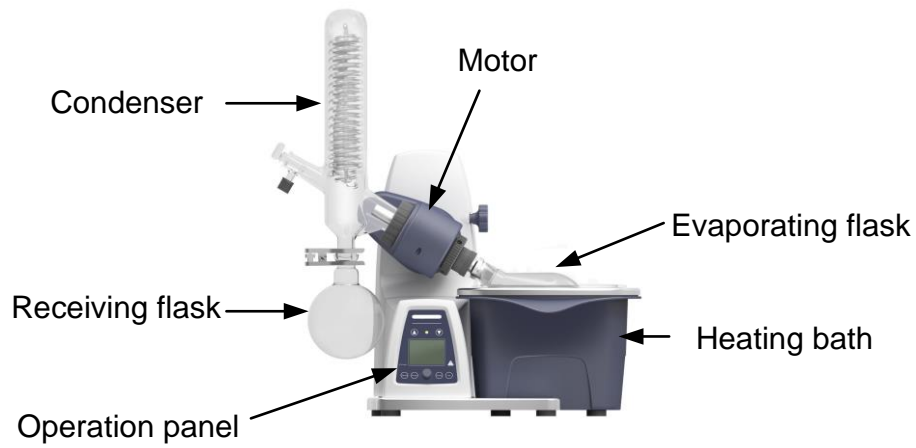
Figure 1

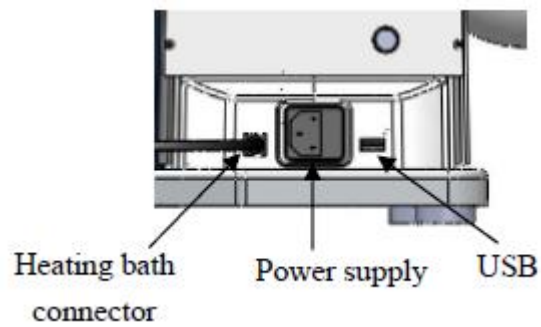
Figure 2

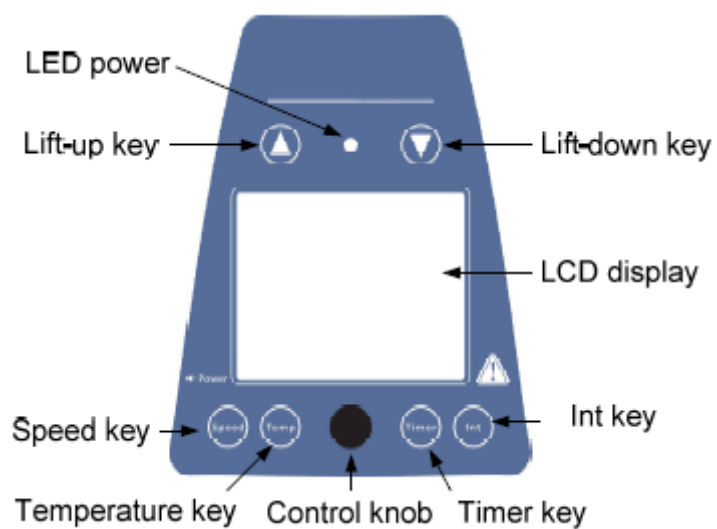
Figure 3

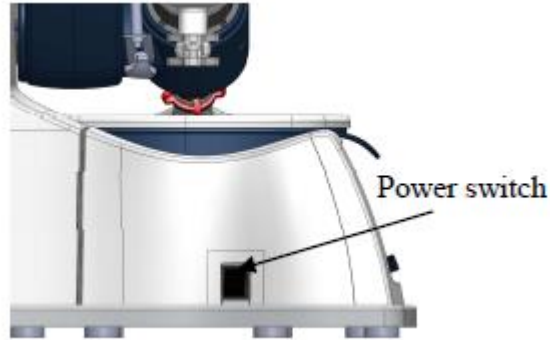
Figure 4

Items	Descriptions
Speed key Speed	Push the key and then rotate the control knob to set the rated speed in the range from 20 to 280 rpm
Temperature key Temp	Push the key and rotate the control knob to set the rated temperature in the range from room temperature to 180°C
Timer key Timer	Push the key and then rotate the control knob to set the rated time in the range from 1 to 999 min
Interval key Int	Push the key and then rotate the control knob to set the interval time for clockwise/counter-clockwise rotation from 1 to 999 s
Control knob	Rotate clockwise to increase the program values. Rotate anti-clockwise to decrease the program values. Press the knob, switch ON/OFF the unit
Lift-up key	Press the key, the lift will go up. Release the key and the lift will stop and remain in the same position
Lift-down key	Press the key, the lift goes down. Release the key the lift will stop and remain in the same position
LCD display	LCD displays the real working state
LED power	Power ON, power LED flashes. When the unit is running, LED power is lit
Power switch	Power ON or OFF

Table 2

4.2 Display

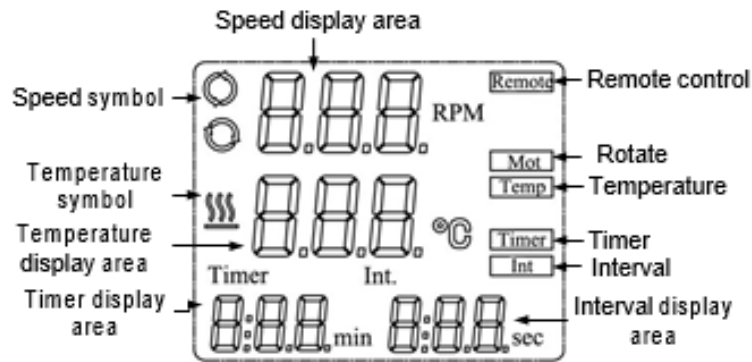
Figure 5

Display	Descriptions
Remote	Remote control
MOT	Rotating function is switched ON
Temp	Heating function is switched ON
Timer	Timer function is switched ON
Int	Interval function is switched ON
Speed display area	While setting the speed, displays the setting value with flashing. The setting value stops flashing when real speed reaches the set point
Speed symbol	Rotating function is switched ON
Temperature display area	While setting the temperature, displays the setting value. When the heating function is switched ON, displays the real value
Temperature symbol	Heating function is switched ON
Timer display area	While setting the timer, displays the rated time. When the timer function is switched ON, it displays the remaining time
Interval display area	Displays the Interval target value

Table 3

5. Installation and operation

This chapter includes how to install the equipment and instructions for a first set-up.

Note: *Inspect the instrument before first use and during unpacking. If necessary, contact the supplier who supplied the equipment and keep the original packaging.*

5.1 Installation

Locate the instrument in a stable place, with a horizontal surface and consider the maximum dimensions of the product. Perform the distillations under vacuum and with the device placed under an extractor hood

5.1.1 Base adjustment

Adjust the base to your workplace by turning the feet as shown in figure 6.

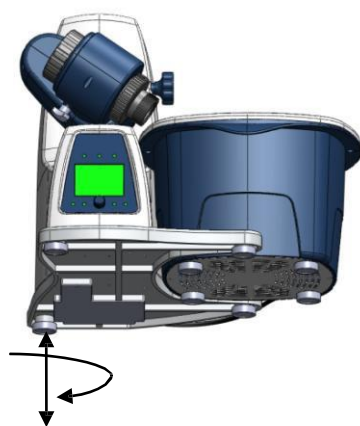
Figure 6

5.1.2 Working position and electrical connections

This equipment has a blocking lock for transport purposes. Proceed as follows so the device adopts its working position:

- Hold the lift with one hand in the same position as the equipment has been provided. Then, with the other hand, unscrew the locking key located on the back of the appliance (A).
- Once the lock has been removed, release the lift and you will see that it begins to rise slowly to the working position. The covered distance is 150 mm.
- Connect the equipment to the power supply (B)
- Connect the thermostatic bath to the device (C)

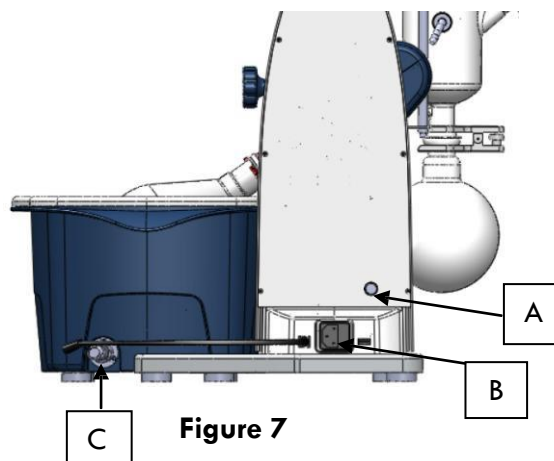
Figure 7

**Attention!**

Make sure the required voltage coincides with electric supply
Connect the instrument to a grounded power source

5.1.3 Adjusting the angle of rotation

- Using the knob (D), you can adjust the rotation angle of the evaporating flask. Turn the key anticlockwise to modify the rotation angle.
- Adjust, by hand, an angle of approximately 30°.
- Once the angle is fixed, close the key (D) by turning it clockwise. In this way, we hold the body that contains the glass parts

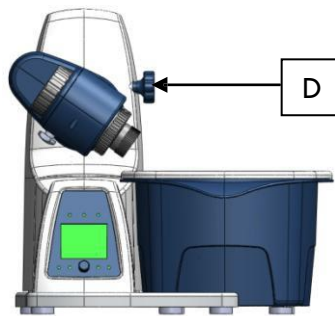
Figure 8

5.1.4 Steam step piece

To insert the evaporator tube (E), proceed with the following instructions:

- Open the locking device (F) by unscrewing the 4 screws located on the main unit by turning counterclockwise
- Insert the evaporator tube until it stops
- Open the locking device on the drive head by turning it 60° counter clockwise (E), (Figure 9)
- Feed the steam pipe in until it stops
- Then lock the locking device by turning it clockwise by 60°
- Close the blocking device by turning it clockwise and inserting again the 4 screws for a better seal

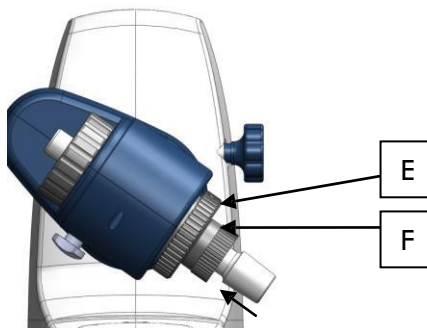
Figure 9

The plastic screw nut (F) helps to create a hermetic seal without tensioning the contact part.

- To install the nut, it is only necessary to insert it through the evaporation tube and screw it in clockwise.

5.1.5 Condenser board

The condenser gasket allows the hermetic seal and guarantees the vacuum of the equipment. It is located in the contact part with the condenser. To install it, observe the following images:

Note: Prior to commissioning, hand-tighten the plastic screw nut (F) left-aligned.

Installing condenser seal (Figure 10 and Figure 11):

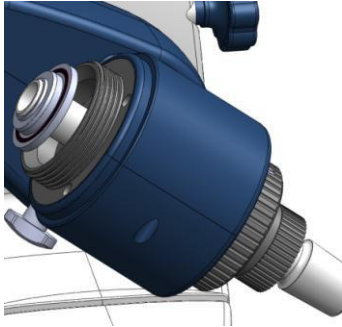
Figure 10

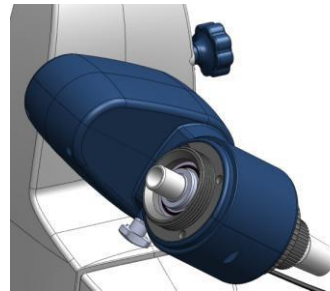
Figure 11

Note: Chapters 5.1.2 and 5.1.3 they do not require installation because they have already been mounted at the factory. It is only necessary to perform this installation when you want to make a change of accessories in the equipment or maintenance. See list of accessories in chapter "9: Accessories and spare parts".

5.1.6 Condenser

- Insert the nut (H) first and place the locking spring (I) into the condenser
- Once the assembly is done, place and install it in the main body by turning the nut (H) clockwise
- To make a better seal, turn on the equipment and rotate at 200 rpm for 2 minutes.

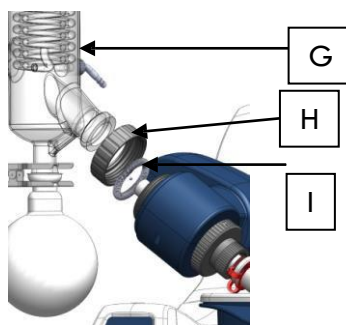
Figure 12

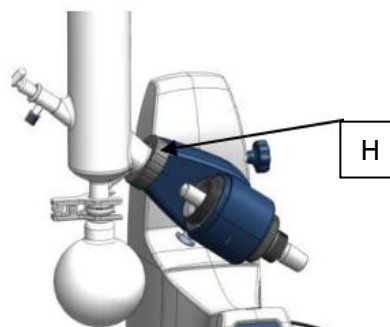
Figure 13

5.1.7 Support rod

The support rod contains a tape that holds the glass assembly. This accessory is optional and has the sole purpose of providing additional stability to the equipment. Especially, we recommend it for glass assemblies that contain a condenser.

-To install the support rod (J), it is necessary to use the locking key (A) removed in chapter "5.1.2 Working position and electrical connections" (Figure 7)

-Position the support rod between the refrigerator and the main body; use the lock key (A) and the screw at the bottom of the motor where there is a slot for that purpose. Once the position is fixed, use the tape provided with all the equipment to surround the condenser and join it with the locking device

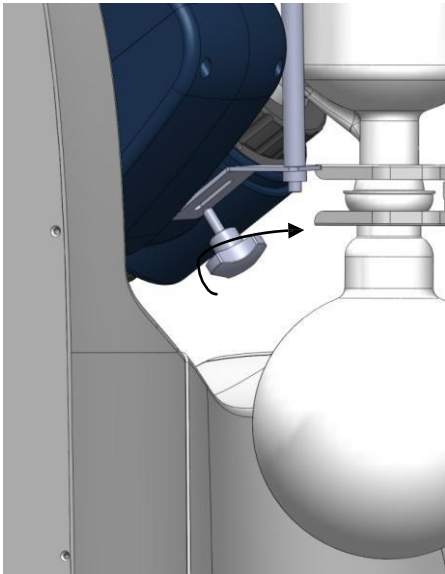
Figure 14

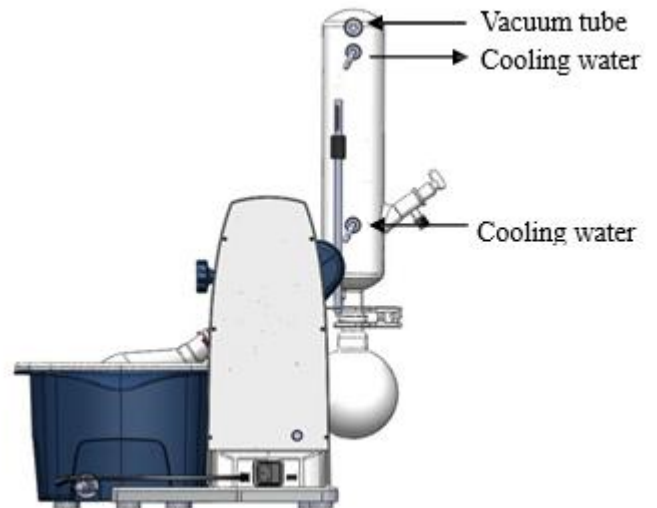
Figure 15

5.1.8 Connections of the tubes

The following image shows a diagram of the connection of the equipment tubes:

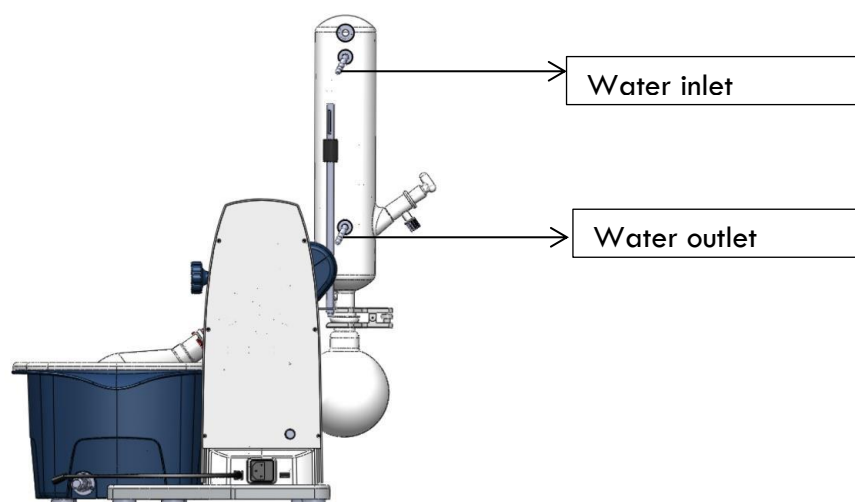
Figure 16

5.1.8.1 Connection pipes with water

While the tubes for condenser water are being connected, keep the following notes in mind:

- Use the connectors for red tubes that we provide with the equipment
- The tubes used must have a diameter equal to the connector, approximately 8mm, see section "9: Accessories and spare parts" to see our references
- For safety reasons, try to secure the tubes with clamps, we recommend that you look at section "9: Accessories and spare parts"
- Check the tubes from time to time and replace them when they are brittle or broken

5.1.8.2 Empty tubes

When connecting the tubes to perform the vacuum in the equipment, keep in mind the following indications:

- Use the connectors for red tubes that we provide with the equipment
- The tubes used must have a diameter equal to the connector, approximately 8 mm, see section "9: Accessories and spare parts" to see our references
- Keep the vacuum tubes as short as possible
- The vacuum pipe does not need to be secured
- Check the tubes from time to time and replace them when they are brittle or broken

5.2 Operation

This chapter explains how to operate the equipment with the different elements and the possible modes of operation properly and safely.

	<p>Warning</p> <p>Risk of injury</p> <p>Do not operate this instrument if the glass is broken</p>
---	---

5.2.1 Motor height adjustment

For a more ergonomic working position, the height of the motor can be adapted to the user requirements. If you want to modify the height, proceed as follows:

- Once the equipment is turned on, press the (up) or (down) button and the height will change
- If you want to finish the height adjustment, stop pressing the button and the movement will stop automatically

5.2.2 Bath temperature adjustment

As soon as you turn on the device with the on / off button, the thermostatic bath will start to heat up if the actual temperature is below the set temperature.

To adjust a temperature, press the "temp" button and, turning the control knob, select the needed temperature. To confirm the action, press the control knob.

	<p>Warning</p> <p>The thermostatic bath can reach temperatures above 100°C. To avoid burns, consider:</p> <ul style="list-style-type: none">- To remove the evaporating flask from the bath- To make sure that the liquid does not overflow the bath when the bottle is submerged
---	---

5.2.3 Adjusting the angle of rotation

The immersion angle inside the thermostatic bath is, by default, 30°. If you need to operate with another angle, for example, when operating with a bottle of different dimension, the angle can be adjusted:

-Turn off the instrument

-With one hand hold the glass assembly, that is, the condenser next to the collecting flask; with the other hand, turn and open the key located on the motor (see chapter 5.1.3).

-Once the key is open, adjust the angle that best suits you. When the angle is adjusted, close the key to fix the rotation position.

	<p>Warning Risk of injury -Never operate this instrument if the glassware is broken</p> <p>Attention! Risk of damage to the instrument -Do not adjust the angle of rotation while the equipment is operating</p>
---	--

5.2.4 Speed adjustment

Perform the following instructions if you want to change the rotation speed of the equipment:

-Press the "speed" button and, by turning the control knob, select the required speed. Press the control knob to confirm the action.

	<p>Warning Risk of damage to the instrument -Adjust the speed of rotation to the dimensions of the bottle and the properties of the solvent</p>
---	---

5.2.5 Timer setting

If you wish to work during specific periods of time, proceed as follows:

-Press the "timer" button and choose the desired operating time (min) by turning the control knob.

-Press the control knob to confirm

-To turn off this function, set the time value to "0" and press the control knob to confirm

5.2.6 Distillate conditions

To achieve optimum distillation conditions, the distillation energy that is provided by the thermostatic bath must be removed by the condenser. To ensure this, operate the instrument according to the following standard rule:

-Cooling water: max. 20°C

-Steam: 40°C

-Thermostatic bath: 60°C

How are these conditions achieved?

- Set the thermostatic bath to 60°C
- Set the cooling temperature to a maximum of 20°C
- With a gauge, grant a speed of approximately 40-50 l/h of cold water through the condenser
- Define the vacuum in such a way that the boiling point of the solvent is 40°C

Note: The corresponding pressure can be found in the table of solvents in chapter 12

Advantages associated with a bath temperature of 60°C:

- The evaporation bottle can be removed without risk of burns
- The rate of water evaporation in the bath is low (low loss of energy)
- The thermostatic bath energy is used with a good degree of efficiency

This rule can also be applied at low bath temperatures, such as:

- Cooling water: 0°C
- Steam: 20°C
- Thermostatic bath: 40°C

5.2.7 Distilled

To start operating with the instrument, comply with the following conditions:

- All electrical connections must be connected
- All sealing is installed correctly
- All joints or contact parts must be greased

To start operating the instrument, follow these indications:

- Turn on the equipment
- Allow cold water to flow through the condenser at a maximum temperature of 20°C and a flow rate of 40-50 l/h
- Set the temperature of the thermostatic bath to the desired value as described above and wait for the bath to reach the target temperature
- Fill the evaporator bottle with the solution to be distilled and make sure it does not exceed a maximum weight of 3 kg
- Assemble the bottle, using tweezers and grease or Vaseline

Note: Choose the pressure so that the boiling point of the solvent is 40°C (see the table of solvents, in chapter 12)

- Adjust the rotation speed

If these operations above are normal, the device is ready to operate. If not, the device may be damaged during transportation, please contact manufacturer/supplier for technical support.

Warning

It is forbidden to transfer the flask while the instrument is in use

6. Faults

This chapter helps to summarize the operations after a minor problem has occurred with the instrument. All possible problems are listed and also, their possible solutions.

6.1 General malfunction and its solutions

The equipment does not turn on:

- Check if the power line is disconnected
- Check that the cable is correctly connected both to the equipment and to the electrical network

The thermostatic bath does not heat:

- Check that it is properly connected to the main body
- Replace the fuse. If the malfunction occurs again, contact the supplier that provided the equipment

The elevator does not work correctly:

- Contact directly with the supplier who provided the equipment

The bottle does not rotate properly:

- Adjust a rotation speed of 0 rpm and, increase the speed of rotation slowly until the rotation motor rotates the bottle
- Restart the instrument

The system is leaking or spilling liquid:

- Grease the joints at the points of contact between materials
- Check the connection of the tubes and adjust them correctly
- The pipe is broken. In this case, replace the tubes
- The sealing system is broken or has not been installed correctly. Check the sealing system explained in chapters 5.1.4 and 5.1.5
- In case you do not fix the fault, replace the vacuum seal or condenser seal

6.2. Malfunction of the vacuum system and its solutions

It does not manage to perform the vacuum in the whole system:

- Please check and make sure that all the joints are greased and perfectly fitted
- If you are using a water vacuum pump to perform the vacuum, open the tap connected to the water vacuum pump to the maximum

The distillation gradually disappears:

- Manually decrease the pressure until the distillation starts again
- If the problem persists, stop the rotation and the pump, open the vacuum valve and check and make sure that there isn't any leaks and that all the joints are greased and perfectly fitted

The distillation stops completely:

- Stop the rotation and the pump, open the vacuum valve and check and make sure that all the joints are greased and perfectly fitted

7. Maintenance and cleaning

This chapter gives you instructions on all maintenance work in order to keep the instrument in good working condition.

Warning

Electrical hazard

The maintenance work must be done with the equipment switched off and remove sources of flammable vapor

7.1 Equipment accommodation

Check the accommodation where you want to place the equipment by default (controls, pests) and clean it regularly with a damp cloth.

7.2 Connection of the tubes

Visually inspect the pipe connection regularly. When the tubes look broken or cracked, replace them with new ones.

“Go to chapter 10 of accessories to see the list of items we can provide.”

7.3 Sealing system

Attention

When removing and reinstalling vacuum seals, ensure not to damage them during handling
Never touch vacuum seals with a sharp object because they may become damaged

7.3.1 Replacement of seals

The seals are subject to a lot of tension and, generally, it is one of the parts that suffer more from the whole team. Check your status regularly and, if necessary, replace them. “See chapter 6.2” to detect possible faults and “chapter 5.1.5” to install the sealing system. Contact the dealer who supplied the equipment for more information.

7.4 Thermostatic bath

The interior surface of the thermostatic bath should be washed if:

- The water used in the bath is calcified or contaminated
- If you use oil instead of water, you must wash the internal surface of the bath when the oil varies in appearance (color, viscosity ...). For this purpose, empty the thermostatic bath in a safe place

7.5 Glass components

To prolong the life of the glass components, consider the following instructions:

-Rinse the glass components with water and a cleaning agent

Note: We advise to wash all glass parts manually.

Note: Use recommended cleaning products in the following situations:

Colorants	Isopropyl alcohol
Construction materials	Water with surfactants / Isopropyl alcohol
Cosmetic products	Water with surfactants / Isopropyl alcohol
Food products	Water with surfactants
Fuels	Water with surfactants

Table 4

8. Specifications

Items	Specifications
Voltage [VAC]	100-240
Frequency [Hz]	50/60
Power [W]	1400
Motor	External rotor brushless motor
Speed Range [rpm]	20-280
Speed Display	LCD
Reversible Direction of Rotation	Yes
Temperature Range [°C]	RT -180
Heat Control Accuracy [°C]	±1
Temperature Display	LCD
Heat Output [W]	1300
Lift	Motor
Stroke [mm]	150
Timer [min]	1-999
Overall Dimensions [D×W×H, mm]	465×457×583
Weight [kg]	15
Permissible Ambient Temperature [°C]	5 - 40
Permissible relative humidity	80%
Protection Class acc. To DIN E N60529	IP20
USB	Yes

Table 5

9. Accessories and spare parts

This chapter includes accessories, sealing items and glassware that can be provided.

Use only the accessories of LABBOX to guarantee the operation of the equipment and ensure a good functioning of the system.

9.1 Glass assembly

Labbox reference	Description
REVA-A01-001	Vertical glass set, includes condenser, distillation flask (269/32) and collector flask (35/20)
REVA-A02-001	Collecting flask (35/20), capacity 1000ml
REVA-A03-001	Piece steam step

Table 6

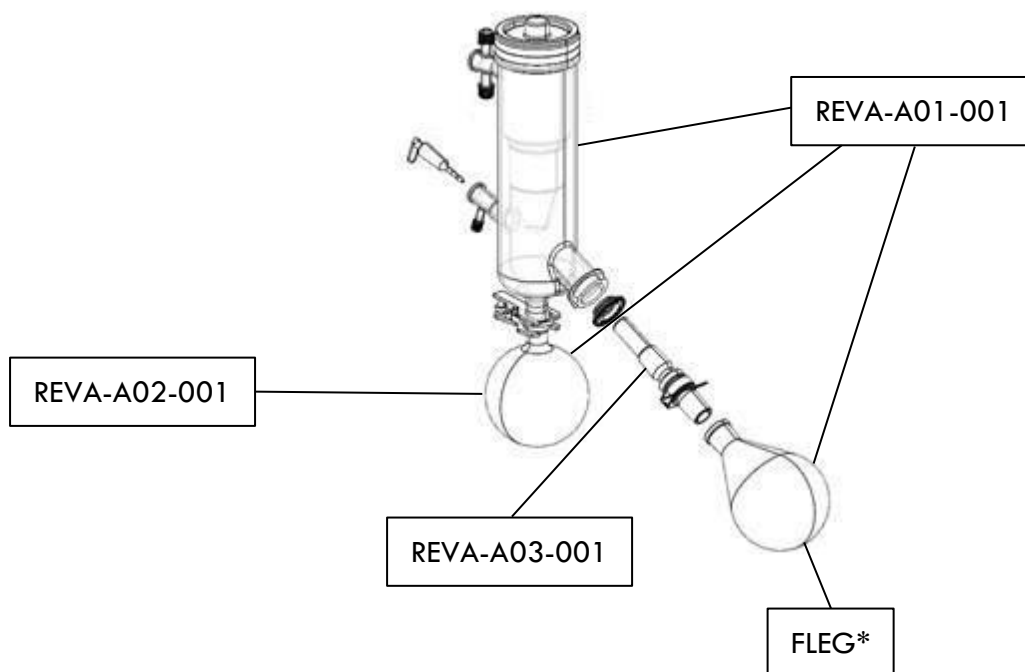
Figure 17

*

Evaporation flask	Capacity
FLEG-051-001	50 ml
FLEG-101-001	100 ml
FLEG-250-001	250 ml
FLEG-500-001	500 ml
FLEG-1K0-001	1000 ml
FLEG-2K0-001	2000 ml

Table 7

9.2 Vacuum references and accessories

Labbox reference	Description
REVA-A04-001	Vacuum seal
REVA-A05-001	Evaporating flask clamp
LTXT-007-005	Latex tubing, amber, for 8/9 hose, 5m
RUTT-008-002	Rubber vacuum tubing for 9/10 hose, 2m
VACU-R20-001	Vacuum pump LBX R20 chemically resistant, 18 l/min, -750 mm Hg

Table 8

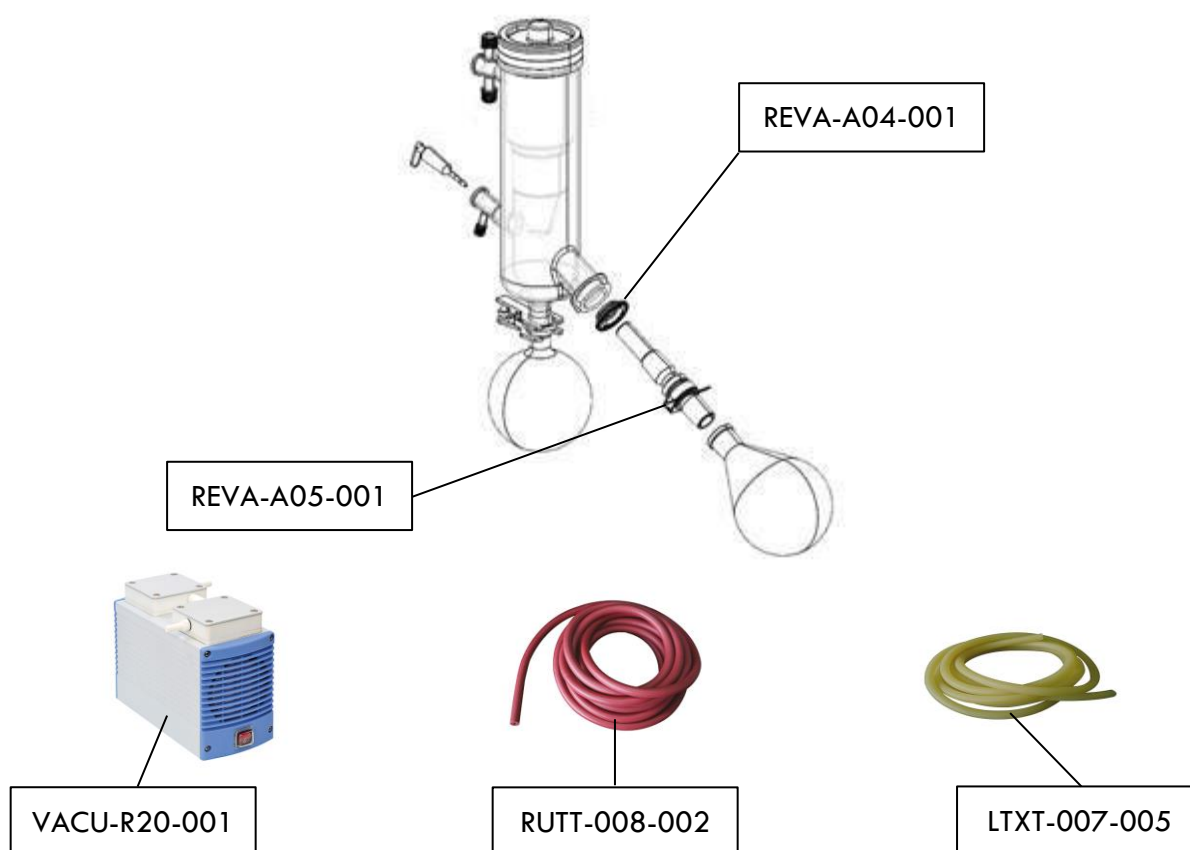
Figure 18

For more information about accessories or laboratory equipment that you need, visit our website www.labbox.com or call +0034 902 99 53 18.

	<p>Attention</p> <ul style="list-style-type: none"> -Electronic devices cannot be cleaned with cleanser -If you require maintenance service, the instrument must be cleaned before shipment to avoid pollution of hazardous substances and also to be sent back in the original packaging. -If the instrument will not be used for a long time, please, switch off and store it in a dry and clean place with a steady room temperature.
---	--

10. Applicable regulations

<p>Manufacture in accordance with the following safety regulations</p>
<p>EN 6110-1 (Safety requirements for electrical equipment for measurement, control and use in laboratories - Part 1 / general requirements) UL 3101-1 (Electrical equipment for laboratory use - Part 1: General requirements) CAN/CSA C22.2 (1010-1) EN 61010-2-10 (Heating)</p>
<p>Manufacture according to the following EMC regulations</p>
<p>EN 61326-1 (Electromagnetic Compatibility)</p>
<p>Associated EU Directives:</p>
<p>EMC Directive: 89/336/EWG Instrument directive: 73/023/EWG</p>

Table 9

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Note: This equipment has been tested to comply with the limits for an electrical device of class A with compliance with part 15 of the FCC rules. These limits are designed to provide protection against interference when the equipment operates in a commercial environment. This equipment generate, uses and can radiate radio frequency energy, therefore, if it is not installed and used in accordance with this instruction manual, it may cause interference in radio communications. Using this equipment in residential areas could cause harmful interferences, in this case, the user will have to correct these interferences on their own.

11. Storage and transportation

This chapter introduces you on how to pack away the instrument, how to prepare it for transport, specific storage and transport conditions.

11.1 Storage and transport

	<p>Alert</p> <ul style="list-style-type: none"> -Remove all hazardous substances from the instrument and clean all parts of the equipment thoroughly -Always keep the equipment in a dry and clean place, with good ventilation and free of corrosive gases -Store and transport the equipment in its original packaging
---	--

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